FUJIFILM**Wako**

Code No. 058-09261 (1mL)

EV-Save™ Extracellular Vesicle Blocking Reagent

EV-Save™ Extracellular Vesicle Blocking Reagent is a polymer reagent to prevent adsorbing extracellular vesicles in cell culture supernatants to labware such as tubes and pipet tips, which reduces the loss of extracellular vesicles during experiments and storage. Add into samples before ultrafiltration of cell culture supernatants, isolation and storage of extracellular vesicles.

The combination of EV-Save™ Extracellular Vesicle Blocking Reagent and MagCapture™ Exosome Isolation Kit PS (Code No. 299-77603, 293-77601) is recommended because of reducing the loss of extracellular vesicles effectively.

【Storage】

Store at -20°C

【Procedure】

〈For isolation of extracellular vesicles from cell culture supernatants using MagCapture™ Exosome Isolation Kit PS (Code No. 299-77603, 293-77601)〉

1. Add EV-Save™ Extracellular Vesicle Blocking Reagent to be diluted 100-fold into cell culture supernatants and mix it by inverting or tapping.
In the case of concentrating cell culture supernatants by an ultrafiltration unit, add EV-Save™ Extracellular Vesicle Blocking Reagent into cell culture supernatants before ultrafiltration.
2. Proceed with an experiment to the step of "Elution of extracellular vesicles" described in the instruction of MagCapture™ Exosome Isolation Kit PS.
3. Add EV-Save™ Extracellular Vesicle Blocking Reagent to be diluted 100-fold into "Exosome Elution Buffer" accompanied with MagCapture™ Exosome Isolation Kit PS and mix it by inverting or tapping.
4. Elute extracellular vesicles according to the instruction of MagCapture™ Exosome Isolation Kit PS.
5. Store the eluted extracellular vesicles, or proceed to the next experiment.

〈For isolation of extracellular vesicles from cell culture supernatants using other methods〉

1. Add EV-Save™ Extracellular Vesicle Blocking Reagent to be diluted 100-fold into cell culture supernatants and mix it by inverting or tapping.
2. Add EV-Save™ Extracellular Vesicle Blocking Reagent to be diluted 100-fold into isolated extracellular vesicles and mix it by inverting or tapping.
3. Store it or proceed to the next experiment.

【Related products】

Code No.	Description	Size
299-77603	MagCapture™ Exosome Isolation Kit PS	2 purifications
293-77601		10 purifications
290-80301	PS Capture™ Exosome Isolation Resin Kit	1 kit (0.5 mL Slurry)
297-79201	PS Capture™ Exosome ELISA Kit (Anti Mouse IgG POD)	96 reactions
297-79701	PS Capture™ Exosome Flow Cytometry Kit	300 reactions
016-27061	Anti CD63, Monoclonal Antibody (3-13)	20 µL
012-27063		100 µL
018-27641	Anti CD63, Monoclonal Antibody (3-13), Fluorescein Conjugated	For 25 tests
014-27643		For 100 tests
290-35591	Magnet Stand	1 each
295-71701	microRNA Extractor™ SP Kit	50 reactions

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
 Telephone : + 81-6-6203-3741
 Facsimile : + 81-6-6201-5964
<http://www.wako-chem.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
 Richmond, VA 23237
 U.S.A.
 Telephone : + 1-804-271-7677
 Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

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FUJIFILM Wako Chemicals Europe GmbH